


FagronLab™ BASIC

Quick - Start


How to register your FagronLab™ BASIC device



- Go to us.fagron.com/en-us/fagronlab.com/re
- Enter the 12 digit serial number
- Enter the four digit unlock PIN on your device

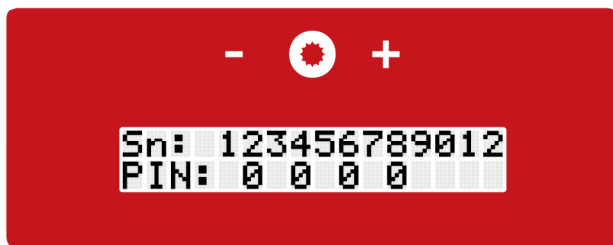


Display handling

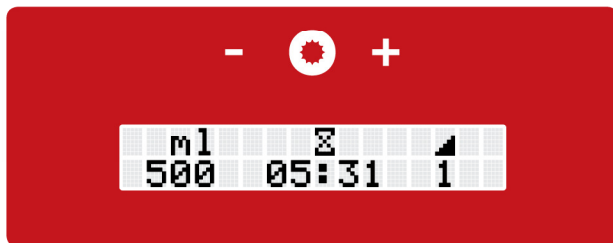

- Touch panel with:
 - Minus button (left)
 - Compounding button (middle)
 - Plus button (right)
- Start display shows jar size, mixing time and mixing speed
- After pushing the compounding button for two seconds "Start" appears on the display and the jar can be mounted

Lift arm with integrated quick-lock coupling

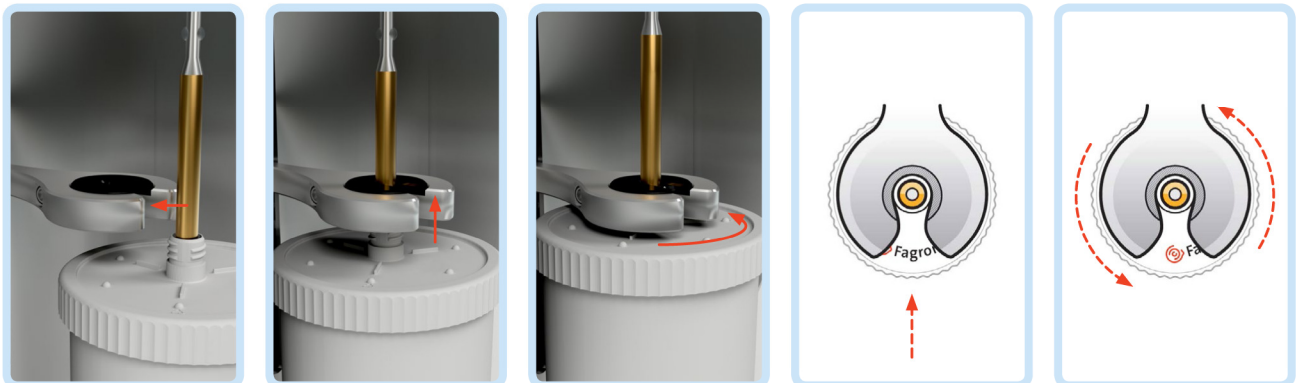

- Hold the FagronLab™ logo on top of the jar to the front; insert the mixing shaft directly from the front into the lift arm
- Insert the mixing unit from below
- Apply slight pressure from the top onto the lift arm. Fix the jar with a slight turn counterclockwise (about a quarter turn) into the lift arm
- Move the mixing unit in the lift arm by hand all the way up and push the compounding button