


FagronLab™ EMP

Quick - Start


How to register your FagronLab™ EMP device



- Go to us.fagron.com/en-us/fagronlab.com/re
- Enter the 12 digit serial number
- Enter the four digit unlock PIN on your device

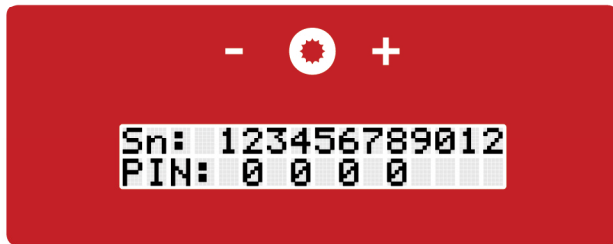

Display handling

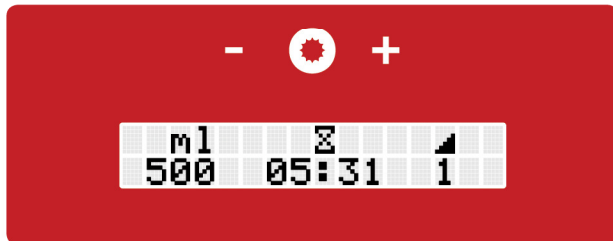

- Touch panel with:
 - Minus button (left)
 - Compounding button (middle)
 - Plus button (right)
- Start display shows jar size, mixing time and mixing speed
- After pushing the compounding button for two seconds the lift arm moves automatically into the mounting position

Lift arm with integrated quick-lock coupling



- Hold the FagronLab™ logo on top of the jar to the front; insert the mixing shaft directly from the front into the lift arm
- Insert the mixing unit from below
- Fix the jar with a slight turn counterclockwise (about a quarter turn) into the lift arm and push the Compounding button