

# FagronLab™ PRO


## Quick - Start

### How to register your FagronLab™ PRO device



- Go to [us.fagron.com/en-us/fagronlab.com/re](https://us.fagron.com/en-us/fagronlab.com/re)
- Enter the 12 digit serial number
- Enter the activation code on your device

## Display handling

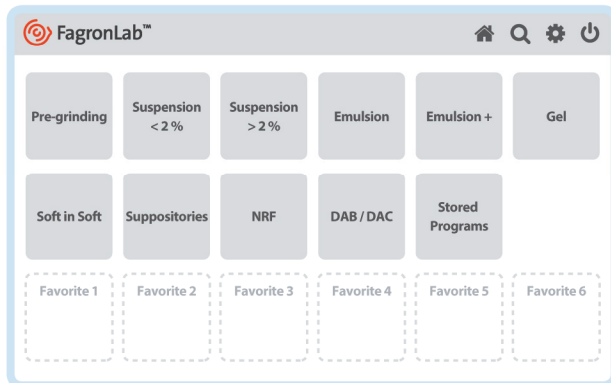

- Touch screen reacting to pressing directly the respective symbols with your finger
- Main display displaying the mixing programs
- After selecting the program, scan of the jar size via the integrated camera or manual input by selecting "jar size"

## Lift arm with quick-lock coupling



- Hold the FagronLab™ logo on top of the jar to the front; insert the mixing shaft directly from the front into the lift arm
- Insert the mixing unit from below
- Fix the jar with a slight turn counterclockwise (about a quarter turn) into the lift arm and push the start button